Auxiliary Display Kit
INSTRUCTION MANUAL

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This Auxiliary Display Kit is for use with OHAUS Scout STX / SPX / SKX / SJX series, Courier 7000 and Defender i-DT25P.

KIT CONTENTS

- Auxiliary Display (PAD7)
- Mounting Brackets (2)
- Instruction Manual
- RS232 Interface
- Power Pack (1) and Plugs (4)

INSTALLATION


1. Set up the Auxiliary Display (PAD7), Power adaptor and RS232 Interface per below.

2. Power on the Auxiliary Display (PAD7).

Install the proper plug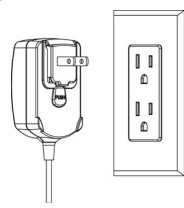



Connect AC Adapter to the proper AC supply

Note: Please use the supplied power adaptor in this kit.

3. Connect the RS232 Interface to the balance.
Install the Interface module on the mini DIN port (round connector) located at the rear side of the balance as shown.

Notes:

- Please make sure the small triangle mark (▲) on the round connector is at the bottom and then plug in.

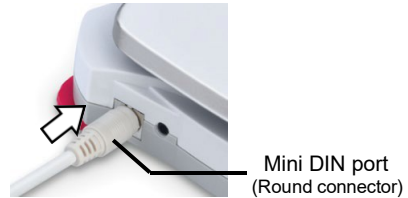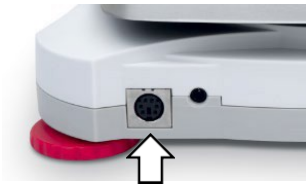4. Turn on the balance and enter the balance's sub menu (Menu/RS232) to configure the parameters as below:

| | |
|------------|--------|
| Baud Rate: | 2400 |
| Parity: | 7 none |
| Stop bit: | 2 |
| Handshake: | none |

5. The Auxiliary Display operates automatically.

If communication is lost, power the balance and the auxiliary display Off and On to restore the connection with the Auxiliary Display.

AUXILIARY DISPLAY MOUNTING & PLACEMENT OPTIONS

- **Table top placement:** The Auxiliary Display may be placed on a flat surface without additional hardware.
- **Direct wall mounting:** The Auxiliary Display may be mounted to a wall using two screws (not included). Select appropriate sized screws that fit into the holes in the back of the Display.
- **Bracket wall mounting/table mounting:** Attach the two mounting brackets. Position the Auxiliary Display where desired on the wall or table.

Note: When installing the mounting brackets, insert round pins into holes in the back of the display and then slide each bracket to lock into place. Mounting brackets may be attached in two different orientations to provide optimum viewing angle.

TECHNICAL DATA

| Model | PAD7 |
|---------------------------|------------------------------------------------------------------------|
| Operating Temperature | 5 °C to 40 °C |
| Relative Humidity | 15% to 80% at 31°C non-condensing, decreasing linearly to 50% at 40°C |
| Altitude | Up to 2000m |
| Storage temperature range | -10 °C to +60 °C |
| Power requirements | 8-14.5V ~ 50/60Hz 40mA or 8-20V $\overline{\text{---}}$ 40mA (min-max) |
| Protection | Pollution degree 2, Overvoltage category II For indoor use only |

Register your accessory under *Warranty Registration* at www.ohaus.com

COMPLIANCE

This accessory has been tested and complies with the approvals listed in the applicable Instruction Manual.

This product conforms to the EMC Directive 2004/108/EC and the Low Voltage Directive 2006/95/EC. The complete Declaration of Conformity is available from Ohaus Corporation.

This device complies with Part 15 of the FCC Rules. Operation is subject to the following two conditions: (1) this device may not cause harmful interference, and (2) this device must accept any interference received, including interference that may cause undesired operation.

This equipment has been tested and found to comply with the limits for a Class B digital device, pursuant to Part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference in a residential installation. This equipment generates uses and can radiate radio frequency energy and, if not installed and used in accordance with the instructions, may cause harmful interference to radio communications. However, there is no guarantee that interference will not occur in a particular installation. If this equipment does cause harmful interference to radio or television reception, which can be determined by turning the equipment off and on, the user is encouraged to try to correct the interference by one or more of the following measures:

- Reorient or relocate the receiving antenna.
- Increase the separation between the equipment and receiver.
- Connect the equipment into an outlet on a circuit different from that to which the receiver is connected.
- Consult the dealer or an experienced radio/TV technician for help.

Please note that changes or modifications not expressly approved by the party responsible for compliance could void the user's authority to operate the equipment.

DISPOSAL

In conformance with the European Directive 2002/96 EC on Waste Electrical and Electronic Equipment (WEEE) this device may not be disposed of in domestic waste. This also applies to countries outside the EU, per their specific requirements.

Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. If you have any questions, please contact the responsible authority or the distributor from which you purchased this device.

Should this device be passed on to other parties (for private or professional use), the content of this regulation must also be related. For disposal instructions in Europe, refer to www.OHAUS.com/weee. Thank you for your contribution to environmental protection.

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P/N 30269077 D © 2016 OHAUS Corporation, all rights reserved / todos los derechos reservados/ tous droits réservés/ Alle Rechte vorbehalten / Tutti i diritti riservati / 版权所有 / 모든 권리 보유 / Все права защищены / すべての権利予約
OHAUS Corporation